



PharmaPure®

Low Spallation Peristaltic Pump Tubing

PharmaPure® Flexibility

PharmaPure® is a premium, low spallation, biologically compatible peristaltic pump tubing developed especially for pharmaceutical, biotechnology, and laboratory applications. This tubing provides excellent pump life with ultra-low particulate spallation and very low gas permeability compared to standard silicone tubing.

Lower Spallation, Lower Risk of Contamination

The superior flex life characteristics of PharmaPure® simplify biopharmaceutical manufacturing processes by reducing production downtime due to pump tubing failures.

Its excellent wear properties allow the product to provide extremely low rates of spallation as compared to silicone and other alternate materials.

Because PharmaPure® has low permeability and lower absorption characteristics, it is ideal for protecting sensitive fluids from gas ingress and concentration changes due to fluid absorption.

Fully Characterized and Biocompatible

PharmaPure® tubing comes complete meeting the requirements of USP <88> Class VI, and/or USP <87>, and/or ISO 10993-5 which can be found in the Validation Guide Summary on the Saint-Gobain Bioprocess Solutions website.

Features / Benefits

- Ultra-low particulate spallation
- Outlasts silicone tubing in peristaltic pumps
- Provides an excellent barrier with very low permeability

Typical Applications

- Cell harvest and media process systems
- Vaccine manufacturing
- Bioreactor process lines
- Aseptic filling
- Diagnostic test products
- Production filtration and fermentation

PharmaPure® Tubing Standard Sizes

Part Number	I.D. inches (mm)	O.D. inches (mm)	Wall Thickness inches (mm)	Length feet (m)	Minimum Bend Radius inches (mm)	Average Burst Pressure at		Max. Suggested Working Pressure at		Vacuum Rating In. Hg (mm Hg)	
						73°F *psi (bar)	180°F *psi (bar)	73°F *psi (bar)	180°F *psi (bar)	73°F (23°C)	180°F (82°C)
AL242606	1/32 (0.8)	5/32 (4)	1/16 (1.6)	25 (7.6)	1/2 (13)	190 (13)	115 (8)	38 (2.6)	23 (1.6)	29.9 (760)	29.9 (760)
AL242002	1/16 (1.6)	1/8 (3.2)	1/32 (0.8)	25 (7.6)	1/2 (13)	100 (7)	65 (4.5)	20 (1.4)	13 (0.9)	29.9 (760)	29.9 (760)
AL242003	1/16 (1.6)	3/16 (4.7)	1/16 (1.6)	25 (7.6)	1/2 (13)	135 (9.5)	90 (6)	27 (1.9)	18 (1.2)	29.9 (760)	29.9 (760)
AL242005	3/32 (2.4)	7/32 (5.5)	1/16 (1.6)	25 (7.6)	1/2 (13)	135 (9.5)	65 (4.5)	27 (1.9)	13 (0.9)	29.9 (760)	29.9 (760)
AL242006	1/8 (3.2)	3/16 (4.8)	1/32 (0.8)	25 (7.6)	3/4 (19)	50 (3.5)	40 (3)	10 (0.7)	8 (0.6)	29.9 (760)	10 (254)
AL242007	1/8 (3.2)	1/4 (6.4)	1/16 (1.6)	25 (7.6)	3/4 (19)	120 (8.5)	60 (4)	24 (1.7)	12 (0.8)	29.9 (760)	29.9 (760)
AL242012	3/16 (4.8)	5/16 (7.9)	1/16 (1.6)	25 (7.6)	3/4 (19)	100 (7)	50 (3.5)	20 (1.4)	10 (0.7)	29.9 (760)	25 (635)
AL242017	1/4 (6.4)	3/8 (9.5)	1/16 (1.6)	25 (7.6)	1-1/4 (32)	75 (5)	35 (2.5)	15 (1.0)	7 (0.5)	29.9 (760)	15 (381)
AL242019	1/4 (6.4)	1/2 (12.7)	1/8 (3.2)	25 (7.6)	1-1/4 (32)	130 (9)	65 (4.5)	26 (1.8)	13 (0.9)	29.9 (760)	29.9 (760)
AL242022	5/16 (7.9)	7/16 (11.1)	1/16 (1.6)	25 (7.6)	1-1/2 (38)	65 (4.5)	35 (2.5)	13 (0.9)	7 (0.5)	29.9 (760)	10 (254)
AL242027	3/8 (9.5)	1/2 (12.7)	1/16 (1.6)	25 (7.6)	1-3/4 (44)	50 (3.5)	30 (2)	10 (0.7)	6 (0.4)	15 (381)	5 (127)
AL242029	3/8 (9.5)	5/8 (15.9)	1/8 (3.2)	25 (7.6)	1-1/2 (38)	95 (6.5)	50 (3.5)	19 (1.3)	10 (0.7)	29.9 (760)	29.9 (760)
AL242038	1/2 (12.7)	3/4 (19.0)	1/8 (3.2)	25 (7.6)	2-1/2 (64)	75 (5)	35 (2.5)	15 (1.0)	7 (0.5)	29.9 (760)	20 (508)
AL242046	5/8 (15.9)	7/8 (22.2)	1/8 (3.2)	25 (7.6)	2-3/4 (70)	60 (4)	30 (2)	12 (0.8)	6 (0.4)	25 (635)	10 (254)
AL242053	3/4 (19.0)	1 (25.4)	1/8 (3.2)	25 (7.6)	3-3/4 (95)	50 (3.5)	20 (1.5)	10 (0.7)	4 (0.3)	15 (381)	5 (127)

*Working pressures are calculated at a 1:5 ratio relative to burst pressure using ASTM D1599.

The values listed for working and burst pressures are derived from tests conducted under controlled laboratory conditions. Many factors will reduce the tubing's ability to withstand pressures including temperature, chemical attack, stress, pulsation and the attachment to fittings. The values shown above are for information only and should not be utilized for specification purposes.


Typical Physical Properties

Property	ASTM Method	Value or Rating
Appearance	--	Opaque White
Durometer Hardness Shore A, 15 Sec	D2240-00	65
Maximum Service Temperature, °F (°C)	--	275 (135)
Low Temperature Embrittlement °F (°C)	D746-98	-89 (-67)
Water Absorption, %, 24 hours @ 73°F (23°C)	D570-98	0.04

Unless otherwise noted, all tests were conducted at room temperature (73°F). Values shown were determined on 0.075" thick extruded strip or 0.075" thick molded ASTM plaques or molded ASTM durometer buttons.



Peristaltic Pump Life Comparison

The table below depicts hours until failure of 1/4" ID x 3/8" OD tubing. In each case, a 4-roller pump head was utilized operating at 300 RPM under room temperature (73°F).

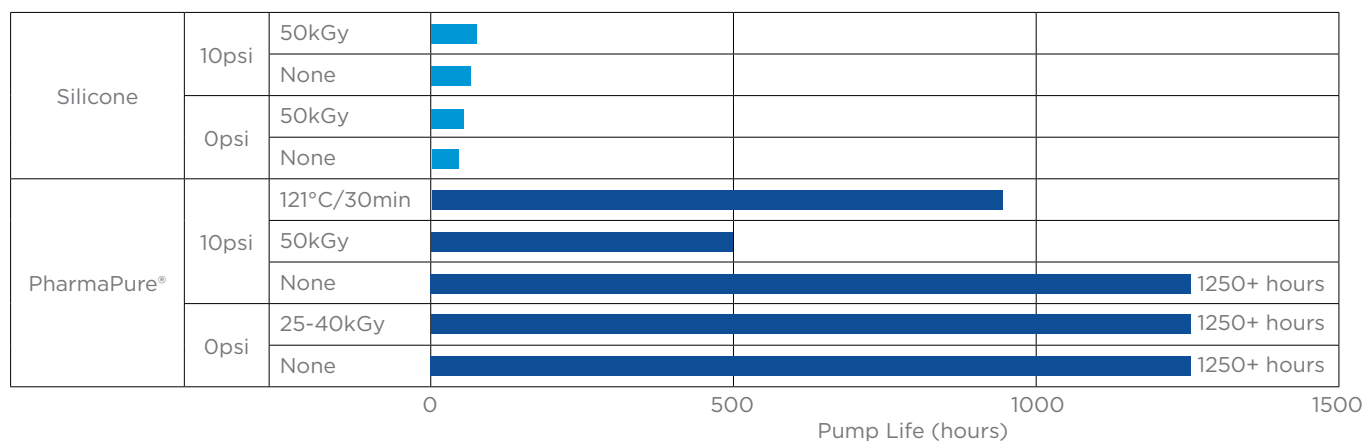

Spallation Comparison

The following test data summarizes the spallation results of select tubing used in a peristaltic pump. In each case 1/16" ID tubing was used in a 2x 6-roller pump head operating at 300 RPM under room temperature (73°F). Spallation is measured with flow imaging cytometry that counts all particles and oil droplets down to 2µm. Results from a minimum of 3 samples were averaged to obtain values

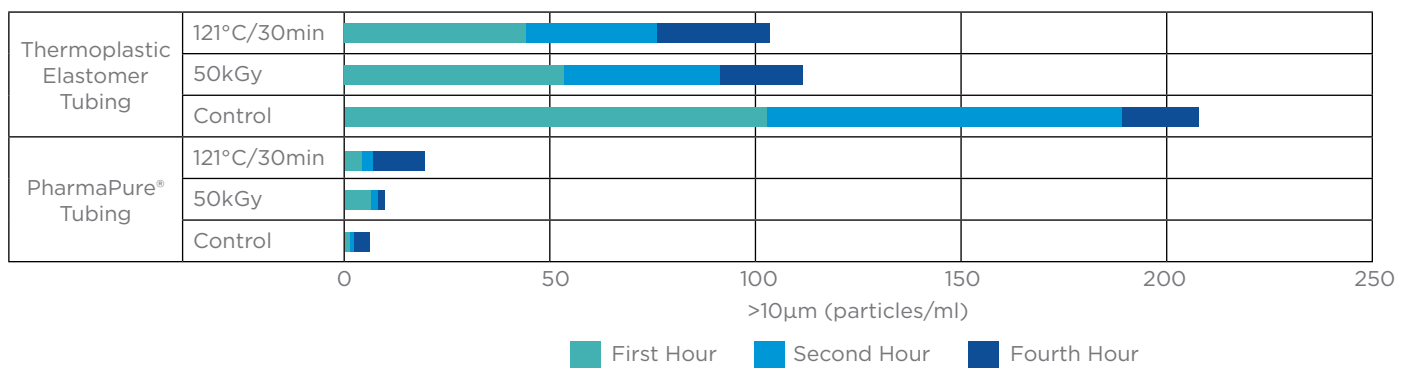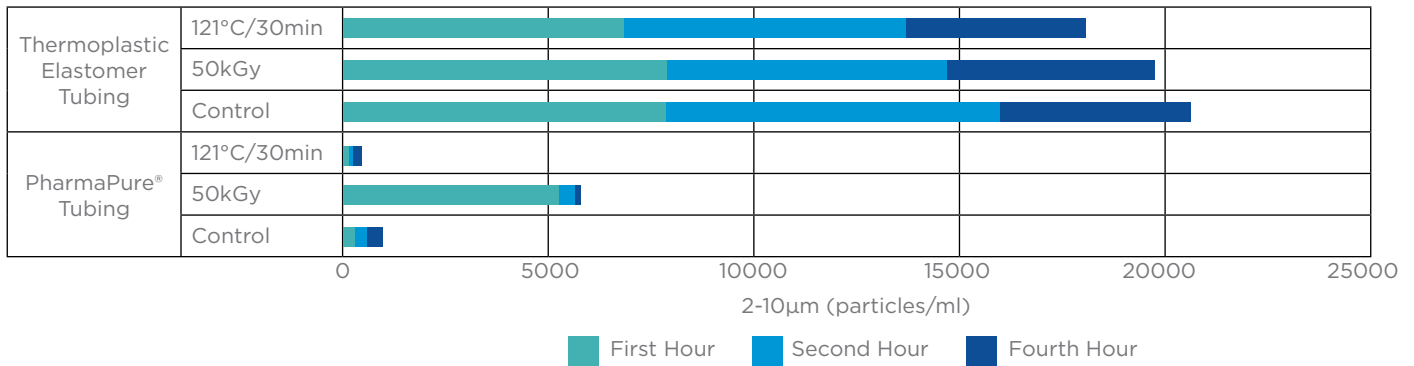

The performance of tubing in peristaltic pumping applications is affected by the conditions of use and equipment utilized, along with size and wall thickness of the tubing tested. The data above is presented for information only and should not be utilized for specification purposes.

Sterilization Methods

Autoclavable	30 min at 121°C
Gas (Ethylene Oxide)	Yes
Irradiation	up to 5.0MRad (50kGy)

PharmaPure® is a registered trademark of Saint-Gobain Life Sciences

Uncontrolled Document - for the controlled version of this document please visit www.biopharm.saint-gobain.com

Contact us today for:

Consultations • Samples • Quotes • Orders



Saint-Gobain Life Sciences

www.biopharm.saint-gobain.com

IMPORTANT: It is the user's responsibility to ensure the suitability, safety and compliance of Saint-Gobain Life Sciences ("SGLS") products/materials for all intended uses and applicable medical regulatory requirements. SGLS assumes no responsibility for product failures due to misuse arising out of the design, fabrication or application of the products into which the materials are incorporated.

WARRANTY: For a period of 12 months from the date of first sale, SGLS warrants this product to be free of defects in materials and workmanship. Our only obligation will be to replace any portion proving defective, or at our option, to refund the purchase price thereof. SGLS DISCLAIMS ALL IMPLIED WARRANTIES OF MERCHANTABILITY AND FITNESS FOR A PARTICULAR PURPOSE.